

## System 75 Telescopic - Aluminium - Wall mounted

Extraction arms	
Wall- and Ceiling mounting 900-1300-3	
Wall- and Ceiling mounting 900-1900-3	
Hoods	
<u>Suction pen 1-7525</u>	
<u>Suction pen 1-7526-5</u>	
Suction nozzle 1-7525-5	75 TE 6
Metal hood 1-7524	
Metal hood 1-7524-5	75 TE 8
<u>Square hood 1-754232-5</u>	75 TE 9
Round hood 1-7528-5	75 TE 10
Round hood 1-7535-5	
Flat hood 1-753324-5	75 TE 12
Brackets	
Wall bracket 2-100-203	75 TE 13
Ceiling columns	75 TE 14
Other accessories	
Reducer 4-100125	75 TE 15
<u>Cover flange 4-250-250</u>	75 TE 16
Support bracket 2-100-5	75 TE 17

# Extraction arm 900-1300-3 – wall- and ceiling





#### **Description**

All components RoHS-compatible

#### Design:


- · Extraction arm for mounting on wall or ceiling
- Working range 950-1300 mm
- Recommended airflow 80 to 180 m<sup>3</sup>/h
- Diameter of tube Ø75/Ø100 mm
- Tubes made of anodized Aluminium surface 10 μ
- Joints made of shatterproof, chemical resistant polypropylene (PP)
- · Colour of joints: white
- · All O-rings maintenance free
- · Integrated valve. When in an open postion out of the flow path to maintain the maximum amount of flow
- All threaded stays, springs and thumbscrews made of acid-proof stainless steel (AISI 316)
- Acid-proof POP® rivets for increased durability in aggressive environments
- · All hoods can be provided with a protective netting (accessory) to reduce the risk of extracting foreign objects
- · Dismantling of the arm without tools for cleaning

Links: Pressure drop chart, Mounting, User's manual, Capture efficiency

### **Working area**

We do not recommend a stationary working position in the upper or lower working area.

All units in mm

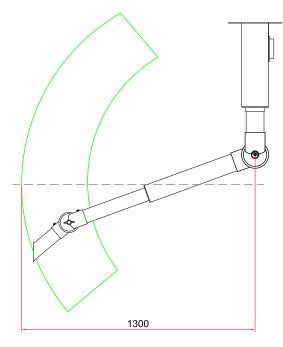



## Extraction arm 900-1900-3 – wall- and ceiling





### **Description**

All components RoHS-compatible.

#### **Design:**


- Extraction arm for mounting on wall or ceiling
- Working range 1300-1900 mm
- Recommended airflow 80 to 180 m<sup>3</sup>/h
- Diameter of tube Ø75/Ø100 mm
- Tubes made of anodized Aluminium surface 10 μ
- · Joints made of shatterproof, chemical resistant polypropylene (PP)
- · Colour of joints: white
- All O-rings maintenance free
- Integrated valve. When in an open postion out of the flow path to maintain the maximum amount of flow
- All threaded stays, springs and thumbscrews made of acid-proof stainless steel (AISI 316)
- Acid-proof POP® rivets for increased durability in aggressive environments
- · All hoods can be provided with a protective netting (accessory) to reduce the risk of extracting foreign objects
- · Dismantling of the arm without tools for cleaning

Links: Pressure drop chart, Mounting, User's manual, Capture efficiency

## **Working area**

We do not recommend a stationary working position in the upper or lower working area.

All units in mm

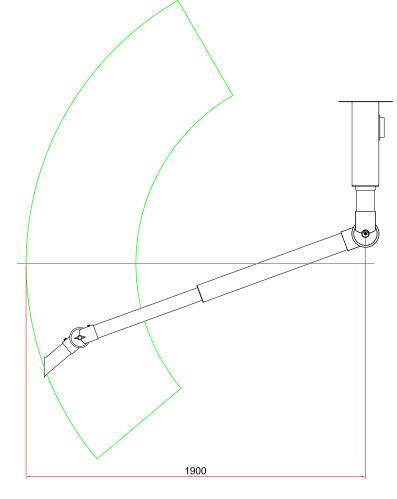








- Suction pen recommended for small concentrated sources of pollution
- High efficiency as the suction pen gets close to the source without obstructing the work process
- Tube made of anodized aluminium surface 10  $\mu$
- Funnel of the suction pen in order to increase the capture efficiency
- · Length: 250 mm



## **Capture efficiency**

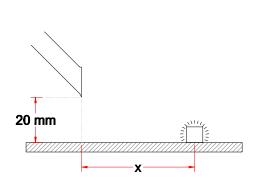

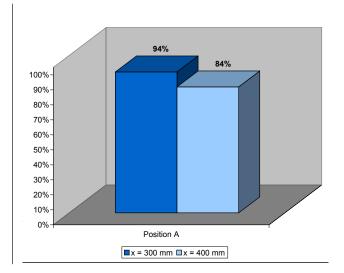

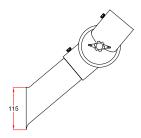

## **Suction pen 1-7526-5**





### **Description**


- Suction pen recommended for small concentrated sources of pollution
- High efficiency as the suction pen gets close to the source without obstructing the work process
- Tube made of anodized aluminium surface 10 µ
- Tip made of shatter proof, chemical resistant polypropylene (PP)
- · Tip colour: white
- Funnel of the suction pen in order to increase the capture efficiency
- · Length: 250 mm



## **Capture efficiency**

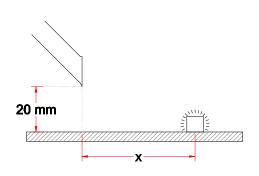

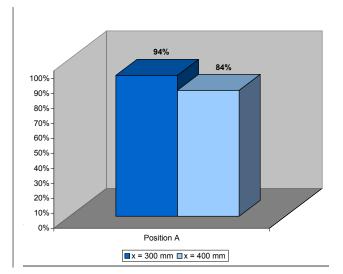

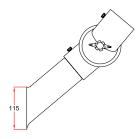

## Suction nozzle 1-7525-5




### **Description**

- Suction nozzle recommended for small concentrated sources of pollution
- Width: 250 mm
- Tube and nozzle made of anodized aluminium surface 10  $\mu$  and shatter proof chemical resistant polypropylene (PP)
- Internal distribution tube in order to increase the efficiency
- · Colour: white



## **Capture efficiency**

Measurements for this product are comparable with article no. 1-5020

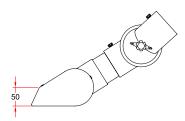





- Metal hood recommended when extracting hot gases, fumes, light dust concentrations etc.
- Suitable for autoclaving
- Diameter of hood: Ø200 mm
- Hood, connection tube and flange made of anodized aluminium surface 10  $\mu$




## Capture efficiency

The relatively low efficiency indicates, that the illustrated position is not optimal for this hood. For system 100 other measurements have been made for a similar round hood - see article no. 1-10024.

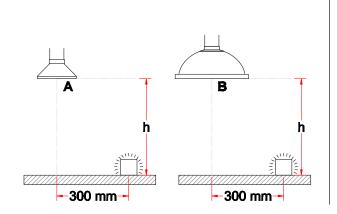

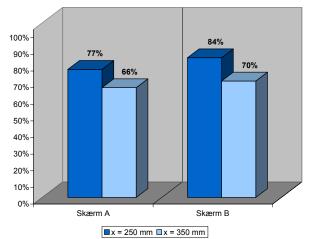

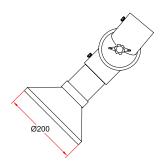





- · Round hood recommended when extracting hot gases, fumes, light dust concentrations etc.
- · Diameter of hood: Ø200 mm
- · Hood made of aluminium with a polyester powder-coating
- Connection tube made of anodized aluminium surface 10  $\mu$
- · Colour of hood and flange: white




## **Capture efficiency**

The relatively low efficiency indicates, that the illustrated position is not optimal for this hood. For system 100 other measurements have been made for a similar round hood - see article no. 1-10024.

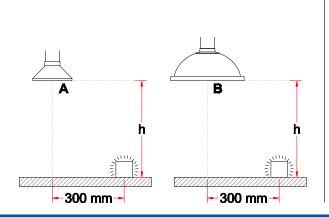

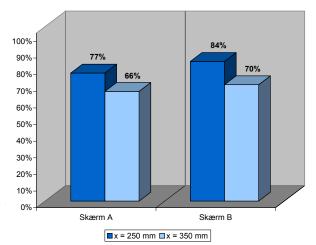

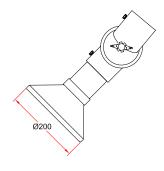

## **Square hood 1-754232-5**





#### **Description**


- Square hood recommended for light and heavy fumes and gases
- Dimension of hood: 420x320 mm
- · Increased stability when moving the hood due to reinforced rim of the hood
- Increased efficiency when placed vertically on a surface
- Hood made of transparent PETG, resistant to solvents
- Connection tube made of anodized aluminium surface 10  $\mu$
- The transparent hood assists in keeping a good view of the work process
- Colours of flange: white



## **Capture efficiency**

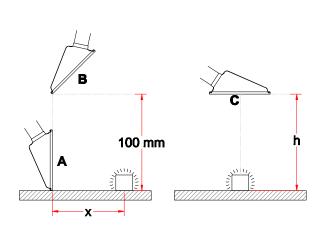

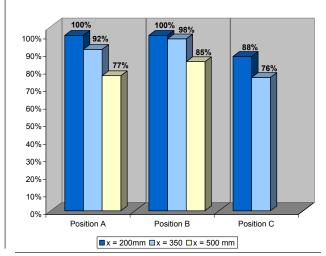

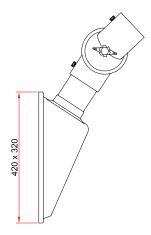



- Round hood recommended for light fumes, gases and over small open vessels
- · Diameter of hood: Ø280 mm
- · Increased stability when moving the hood due to reinforced rim of the hood
- · Increased efficiency at an angled position
- Hood made of transparent PETG, resistant to solvents
- Connection tube made of anodized aluminium surface 10  $\mu$
- The transparent hood assists in keeping a good view of the work place
- Colour of flange: white



### **Capture efficiency**

Measurements of this product are comparable with article no. 1-6328.





- Round hood recommended for light fumes, gases and over small open vessels
- · Diameter of hood: Ø385 mm
- · Increased stability when moving the hood due to reinforced rim of the hood
- · Increased efficiency at an angled position
- Hood made of transparent PETG, resistant to solvents
- Connection tube made of anodized aluminium surface 10  $\mu$
- The transparent hood assists in keeping a good view of the work place
- · Colour of flange: white




### **Capture efficiency**

The relatively low efficiency indicates, that the illustrated position is not optimal for this hood. For system 100 other measurements have been made for a similar round hood - see article no. 1-10050.

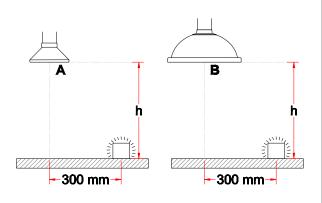


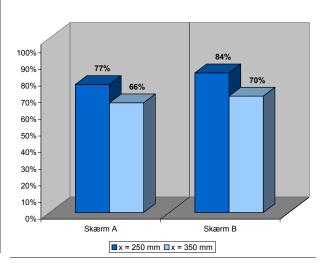

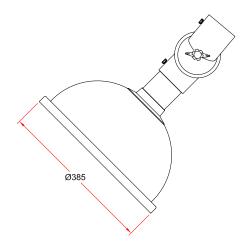






- Flat hood recommended for heavy gases and fumes
- Dimension of the hood: 330x240 mm
- Increased efficiency when placed vertically on a surface
- · Hood made of transparent PETG, resistant to solvents
- Connection tube made of anodized aluminium surface 10  $\mu$
- The transparent hood assists in keeping a good view of the work place
- Colour of flange: white

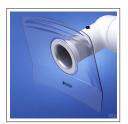

## **Capture efficiency**

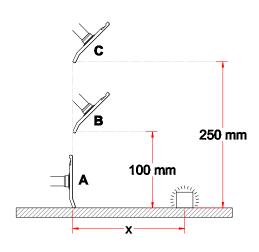

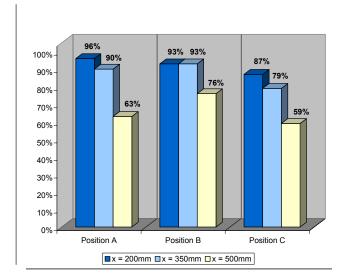


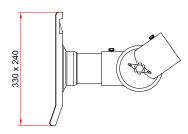




- Wall bracket to mount extraction arms for wall or ceiling mounting to the wall
- Made of steel with a polyester powder-coating on all surfaces or acid proof stainless steel (AISI 316L)
- Colour: white
- Reducer to duct Ø125mm included



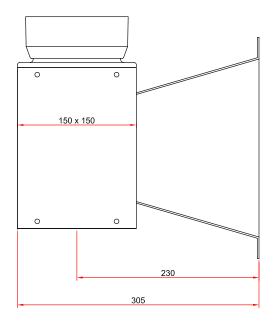

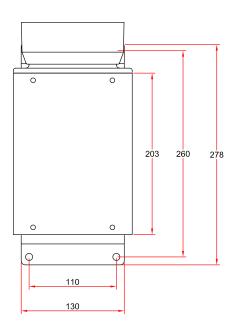



## **Ceiling columns**



## **Description**

- Ceiling column to mount extraction arms for ceiling mounting to the ceiling
- Strong construction made of steel
- Increased durability due to polyester powder-coating on all surfaces
- Side connection: Ø125mm
- · Available with top connection on request
- · Colour: white
- Dimension: 150x150 mm • Length: 500 - 2500 mm

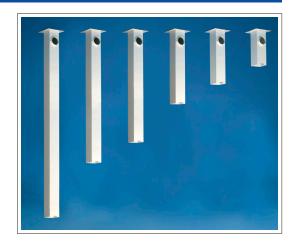


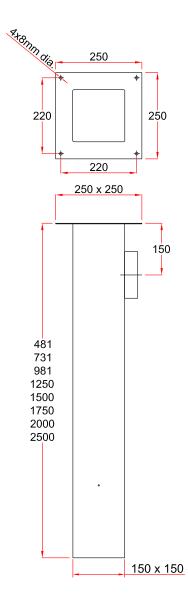



- Reducer to connect extraction arm and ventilating duct
- Reduces from Ø125 Ø100
- Made of polypropylene (PP)
- Colour: White



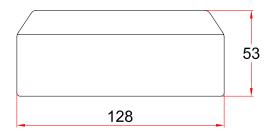

## **Cover flange 4-250-250**



## **Description**

- · Cover flange to mount on false ceiling. Hides the carrying of the ceiling column through the false ceiling
- Dimensions: 250 x 250 mm
- Material: polystyrene
- Fasten with glue or small screws
- · Colour: white



#### **Drawing**

For this product there is no drawing


## **Support bracket 2-100-5**



## **Description**

- For extra support of long ceiling columns when mounted to ceiling racks or with wires to the
- Made of steel with a polyester powder-coating on all surfaces
- · Colour: white
- Dimensions: 250 x 250 mm



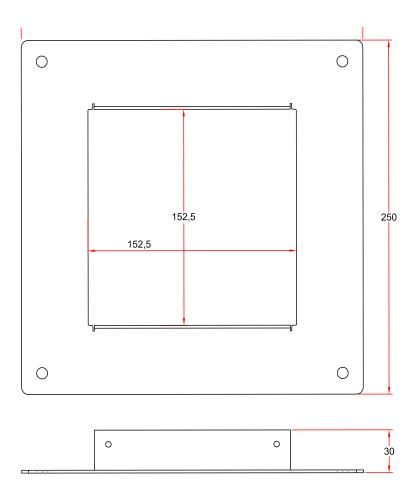

# **Protective netting 5-17**



## **Description**

- Protective netting to be placed in the hood
- Protects against extraction of foreign objects
- Polypropylene (PP)
- Colour available: white



